



English


Deutsch

Français

Italiano

Contact information:
email: spycamera.expert@hotmail.com
Skype: [nastya.huang](#)
Phone: +86 186 6434 1559

Wifi Box Digital Microscope (user manual)
Customer services:spycamera.expert@hotmail.com

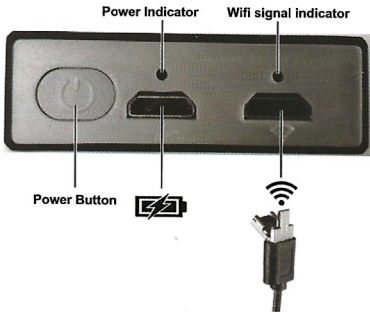


How to charge the wifi box:

Connect the charging port with the socket charger or PC usb port. When the Battery LED stays on, it is charging. When the Battery LED goes off, it is full charged.

1.iPhone/iPad/Android mobile installation (Use wifi box)

Search “wifimage” on App store/Google Play or other app markets. Or you can scan these codes to get the downloading page:




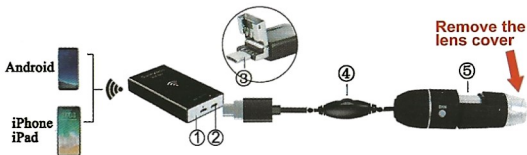

Long press the ①Power button of Wifi box to activate it. Battery red led lights up first. After 5s, WiFi Signal green led lights up. (if not, please charge the wifi box firstly) Connect the microscope's ③micro USB port to the ②Box's WiFi data transfer port. Microscope led will light up. (if not, please check the cable led switch of the microscope).

Go to mobile's WiFi listing and find the WiFi whose name is "coolingtech-xxxx". Then go to the app "wifiimage" and find the device list, rotate the ④LED Control to adjust the light and ⑤Focus Wheel to focus. (must remove the lens cover, or the image can't be clearly)



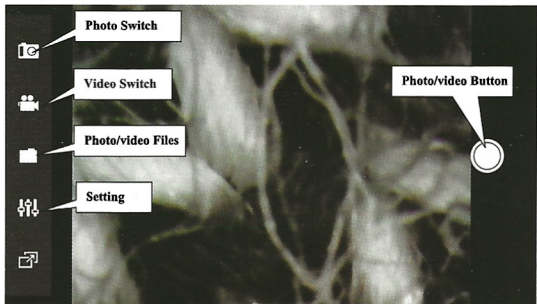

Note:


- a) Low power may cause video transfer interrupted. Please full charge it before using.
- b) If you only see black screen on "wifimage" app, please go to the settings to allow "wifimage" app to use Wifi Data. And the "coolingtech-xxxx" wifi should keep connecting during using.
- c) **SNAP** and **ZOOM** button can only use in **Windows Amcap driver**. In other devices, please use the **focus wheel** to zoom and software's screen-saving function to snap.

2.PC installaiton (Wifi box is unnecessary on Windows or system below) windows 8/10 systems


Step1: Connect the microscope like picture



Step 2: Click Camera

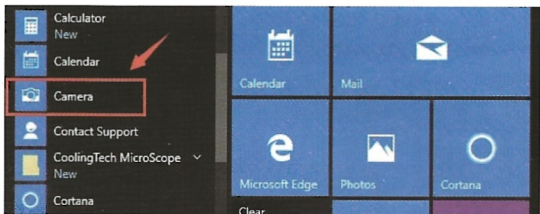

Picture 1

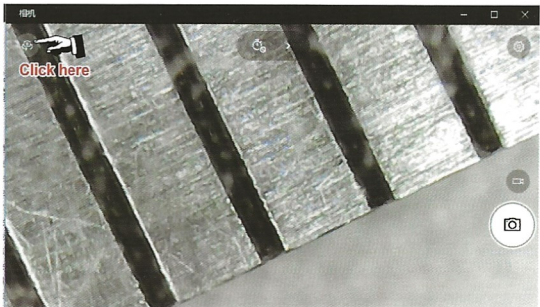


Picture 2


Step 3: Successfully! use the focus wheel(5) to focus

3.How to install the Amcap software in PC / Windows 7/8/10 system(only use this software the ZOOM and SNAP button can work normal,and use this method you needn't run the CD)


Connect the USB port to the PC and then click this link: www.gloptics.net

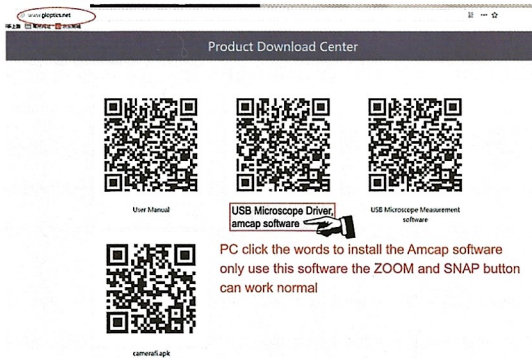

The screenshot shows the 'Product Download Center' page. At the top left is the 'www.gloptics.net' logo. Below it are navigation icons. The main content area features five QR codes arranged in two rows. The first row contains three QR codes with labels: 'User Manual', 'USB Microscope Driver, amcap software' (with a hand icon pointing to the text), and 'USB Microscope Measurement software'. The second row contains one QR code labeled 'camera4apk'.

Click"driver"---and double click"setup.exe"----"Next"--"Install"---"Yes"---"Finish"After you installed the Amcap software,this item  can be appeared on the desktop.Click it, and can be use the microscope normal.

If you want to save the photos when observing,pls click the“Capture”

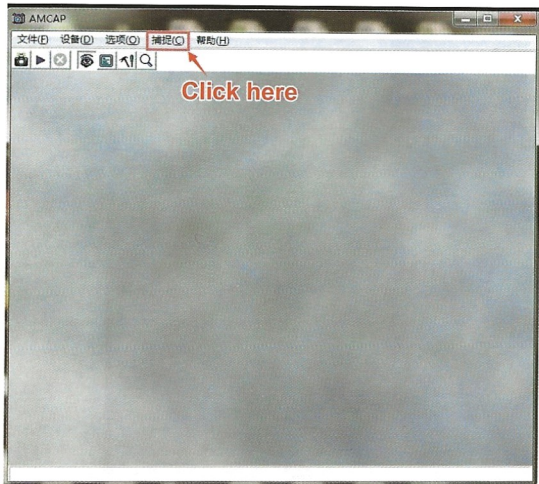

Then click “Storage location of photos”,select the location that you want to save,double click the “Snap”button on the microscope,and then you will find the photos in the place where you just set.

4. How to use on MAC

Mac user can directly use the built-in webcam



5. How to use on Chrome OS

Sunnywoo microscopes can be used with the default Camera app that comes with Chrome. Some Chrome machines with built-in webcam cameras may need to have the device pointer changed. This can be done by going to **Settings > Advanced Settings > Privacy > Content Settings > Camera**, and change the device in the dropdown to "USB Microscope"

Frequently asked questions:

1. Q: Can't find WIFI hot spots?

A: Please fully charge the WIFI box first, then press the power button for 5 seconds and then the green light will appear. Open the mobile WIFI search for "coolingtech XXXX"

2. Q: Can't connect to WIFP?

A: Users of Android system must first turn off "WLAN Assistant" in the WIFI setting option of the mobile phone. If the "WLAN Assistant" option is not turned off, the WIFI hot spot can not be connected .

3. Q: What to do with the problem that the lens is hard to focus?

A: It is recommended to use the PC Amcap software or the mobile phone App: CameraFi (you can download on the site: www.gloptics.net, just scan the "camerafi.apk" code)

4. Q: What to do if the software running gets slower or not smooth during using?

A: It is recommended to restart the computer or uninstall and re-install the amcap and measuring tool.

5. Q: Why my iphone or ipad/Android can't see the image/black screen?

How can I save images/videos on ipad/iphone You need to allow the "wifimage" to use the data. or you can go to settings to set up. IOS user can set up via "Settings-wifimage-Allow on Photos/WLAN&Cellular Data"

6. Q: Why I can't use the ZOOM and SNAP button on mobile phone?

The ZOOM and SNAP button can be use on PC, but not mobile phone, and ZOOM and SNAP also can't be use in measurement software.

If you have any questions, please contact us before write a review, we will provide 100% satisfactory solution.our

Email is: spycamera.expert@hotmail.com

Skype: nastya.hunag